

VACUU·SELECT®

NEW VACUUM CONTROLLER FOR CHEMISTRY AND LIFE SCIENCE

time-saving, easy-to-use, versatile

Asynt



Technology for Vacuum Systems

SIMPLY UNDER CONTROL

VACUU·SELECT® - THE NEW VACUUM CONTROLLER

Looking for an intuitive, easy-to-use vacuum controller? VACUUBRAND has introduced the VACUU·SELECT controller - an application-based controller with state-of-the-art touch screen display for all your vacuum processes.

VACUU·SELECT has you covered. Use manual set-point-control for simple processes, run fully automated distillations or create your own application with simple drag-and-drop editing. VACUU·SELECT comes with pre-defined applications for all common tasks in the lab.

The controller takes care of all vacuum processes for you. For solvent evaporation, VACUU·SELECT detects boiling points and adjusts the vacuum process as necessary. You can virtually forget about the vacuum process and devote your attention to other tasks.

VACUU·SELECT - ready for your process.





SIMPLY EFFICIENT

- intuitive user interface
- pre-defined procedures for individual applications
- easy drag-and-drop editing for creating new procedures
- fully automatic evaporations at the touch of a button
- integrated contextual help
- choose from 14 languages



DESIGNED FOR THE CHEMISTRY LAB

- for all chemical and life science processes
- all wetted parts made from chemical resistant materials
- touch screen display can be operated even with safety gloves
- installation flexibility: put the control panel where you need it
- ready for modern laboratory networks




A SOLUTION THAT ALWAYS WORKS

- as part of VACUUBRAND chemistry pumping units
- independent controller for existing vacuum sources
- fine vacuum control down to 10^{-3} mbar, e.g. for Schlenk lines or lyophilization
- built-in unit for laboratory furniture, vacuum networks and industrial process systems



SIMPLY EFFICIENT

Intuitive user interface

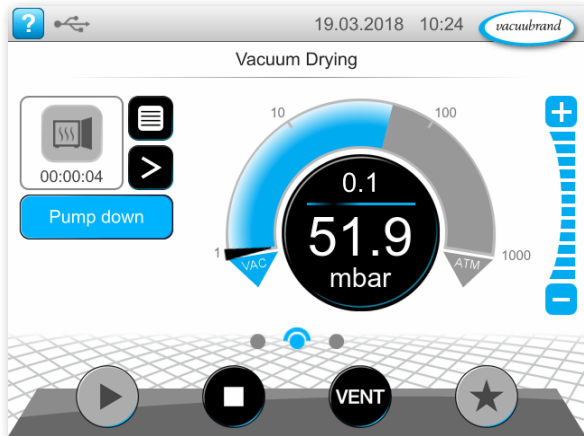

Using the VACUU-SELECT controller is just like using a smartphone. Interactive menus designed specifically for lab applications get you up and running in no time.

application-based navigation

integrated contextual help function with application-specific tips

choose from 14 languages



Ready to use vacuum procedures for all common applications




VACUU-SELECT includes predefined vacuum procedures for all common lab applications. The process steps and control parameters can easily be adjusted to suit any application.

It's as simple as

pick your application

use the predefined parameters or adjust as necessary

press Start

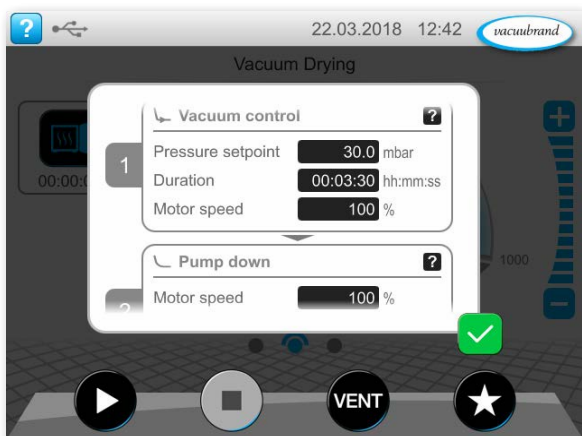



Fully automatic evaporations at the touch of a button




Automatic evaporation mode is perfect for controlling applications like rotary evaporation or vacuum distillation. Simply press the start button and the boiling pressure will be detected automatically and continuously adapted throughout the process. This makes the process run quickly, and without foaming. There's no need for manual adjustments or constant babysitting.

The function "automatic evaporation" works in combination with a **VARIO chemistry diaphragm pump**.



Easy customization through drag and drop editing

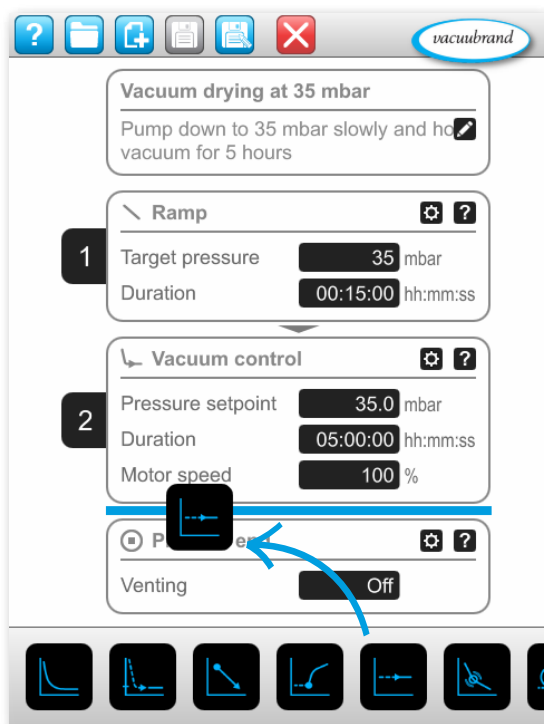

Use the new drag and drop editor to adapt the pre-defined processes for your work. Insert, delete, and modify process steps in an instant to create individualized vacuum processes. Once created, your process will run exactly the same way, time after time. This ensures reproducibility and frees up time for other tasks.

Here's how you create your very own process

drag & drop the desired process steps into the right sequence

adjust the parameters, like target pressure or duration

save your customized application and make it a Favorite

select it from the Favorites menu and hit Start



DESIGNED FOR THE CHEMISTRY LAB

Covers all common vacuum processes



One for all

VACUU-SELECT works with a wide range of vacuum sources - whether integrated into a VARIO pump, used with an existing lab pump, or in connection with a lab vacuum system. All common lab applications are supported by a single controller. VACUU-SELECT simply has you covered.

rotary evaporation, with proven VACUUBRAND fully automatic mode at the touch of a button

all other applications in the rough vacuum range down to 1 mbar

Fine vacuum control down to 10^{-3} mbar, e. g. for Schlenk lines or freeze-drying

vacuum control for kilolabs and pilot plant scale

Durable & chemical resistant



Thick-skinned where it's needed

Proven chemical resistant wetted materials and sealed display electronics protect the controller against corrosion and ensure durability.

glass touch panel display

chemical resistant wetted materials

ceramic vacuum sensor

Touch screen display can be operated even with gloves



You can leave your glove on

Even when wearing thick safety gloves, VACUU-SELECT still responds to your touch.

Variable set up of the control panel



Put the controller where you want it to be

The VACUU-SELECT controller allows you to separate the control panel from the sensor. Leave the pump in the base cabinet and put the display on the bench. Installation flexibility allows you to set up your work space as you like it.

Ready for the future



Your data - always under control

VACUU-SELECT can be integrated into modern networks and data management systems. Read out data using a USB stick or network interface. Modern software allows for simple updates and add-ons.



A SOLUTION THAT ALWAYS WORKS

VACUU-SELECT Complete Controller - for existing vacuum pumps



All-in-one, when and where you need it

The VACUU-SELECT Complete Controller includes the control module, and integrated vacuum sensor, control valve, and venting valve. Available with a table top stand, for lab scaffold mounting, or for panel mounting.

install the controller in your vacuum line, between the pump and your application

the VACUU-SELECT controller detects solvent boiling then holds the pressure level to maintain process control

venting valve allows connection of inert gas

integrated check valve minimizes interference and maximizes vacuum stability



Fine vacuum control configurations down to 10^{-3} mbar



Convenient control down to 10^{-3} mbar

Complete packages provide everything you need for controlling vacuum down to 10^{-3} mbar.

VACUU-SELECT covers fine vacuum applications like freeze drying and Schlenk lines

control range from atmospheric pressure down to 10^{-3} mbar

combination of vacuum controller VACUU-SELECT with fine vacuum sensor, high flow vacuum control valve and all required connections

All-in-one chemistry pumping units

Something great just got better

VACUUBRAND chemistry pumping units - now with VACUU·SELECT



VACUUBRAND chemistry pumping units with valve control

these units combine VACUU·SELECT with VACUUBRAND's powerful, reliable chemistry diaphragm pumps and electronic control valves

the VACUU·SELECT detects solvent boiling then holds the pressure level to maintain process control

independent control for up to two separate processes using a single controller



Speed controlled VARIO chemistry pumping units: unparalleled precision in vacuum control

the ultimate combination of speed controlled pumping and intuitive application-based control minimizes process times and prevents bumping and foaming

pump speed adjusts dynamically to process conditions - running whisper-quiet, minimizing power consumption and service requirements

Built-in vacuum control for lab furniture and vacuum systems



Easy-to-use control for lab vacuum networks

The VACUU·SELECT Complete Controller can easily be built into lab furniture and casework and integrated into VACUU·LAN lab vacuum networks, for convenient, modular, and controlled lab vacuum networks.



ELECTRONIC VACUUM CONTROL

Benefits from electronic vacuum control

- minimizes process times
- enables automated processes
- generates reproducible results
- minimizes risk of foaming and sample loss


State of the art control can make all the difference

VACUU-SELECT supports **two vacuum control procedures**. Classic 2-point control uses a solenoid valve with the VACUU-SELECT controller. Advanced VARIO control achieves unparalleled precision with motor speed control.

Classic 2-point vacuum valve control

Compatible with any vacuum source

- detects the onset of solvent evaporation then maintains a stable vacuum level
- pump is permanently running at full speed

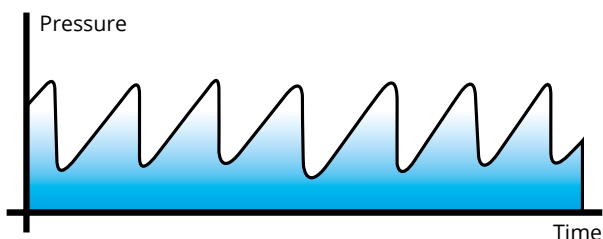


The vacuum controller switches a vacuum valve on and off, holding the vacuum level stable within a tight range or hysteresis. VACUU-SELECT automatically adjusts the hysteresis based on process demands in order to minimize pressure fluctuations and maintain process control. This is especially important for evaporative processes, where the hysteresis is automatically set based on the first vapor pressure detected.

Precise VARIO® control via motor speed control

Available only with VARIO vacuum pumps

- automatic vapor pressure detection and continuous pressure optimization
- matches pumping speed to actual demand at all times

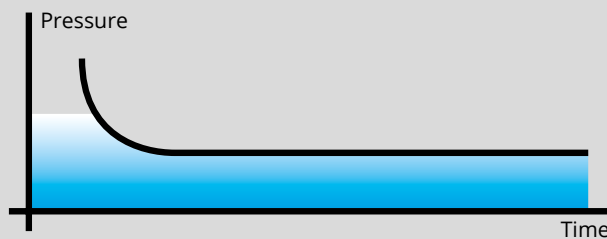

VARIO pumps and pumping units combined with the VACUU-SELECT controller deliver unparalleled precision, free of hysteresis. Pumping speed is continuously and automatically adapted to achieve optimal gas flow based on the rate of solvent evaporation or gas load, and the system volume. VARIO control always seeks to optimize process conditions and minimize process time.

VARIO®

at a glance

Automated evaporation


In Automated Evaporation mode, the new VACUU·SELECT controller automatically detects solvent vapor pressures. With continuous optimization of the vacuum level based on real-time process conditions, there is no need to actively monitor your process - no more babysitting is required. Just press start and walk away.

minimizes process time

reliably maintains control for you - no more babysitting

makes manual adjustments and programming a thing of the past

minimizes foaming and sample loss

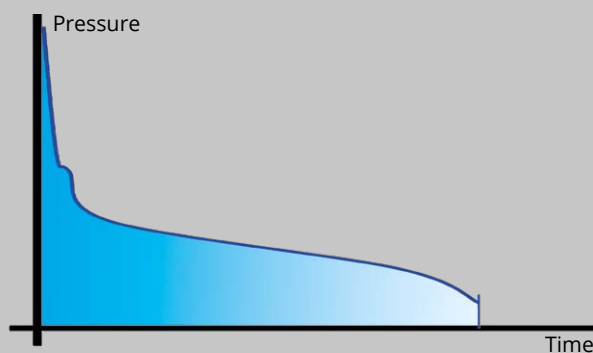

Pump speed on demand

whisper quiet operation

drastically reduced power consumption

decreased service demand

Precisely optimized vacuum

minimizes process times, saving as much as 30% compared to other control methods

reproducible processes means reproducible results



VACUU-SELECT Complete Controller

The VACUU-SELECT Complete Controller is ready for use with existing vacuum sources like pumps or vacuum networks.

The intuitive touchscreen display offers predefined tools for all common vacuum processes. In evaporative processes, the first vapor pressure is detected then the controller uses 2-point control to maintain a stable vacuum. With several other applications included, and the ability to adjust or customize processes to your needs, VACUU-SELECT has you covered for a wide range of lab vacuum applications.

Available in a variety of configurations, VACUU-SELECT integrates seamlessly into virtually any lab. The VACUU-SELECT Complete Controller includes an integrated ceramic vacuum sensor, a vent valve, vacuum control valve, and check valve. This compact unit is easy to connect to your vacuum source.



control range	measuring principle	vacuum connector	Ord. No.	
atm. - 0.1 mbar	capacitive measurement via ceramic diaphragm	hose nozzle DN 6/10 mm	table top:	20700070
			scaffold mounting:	20700080
			built in:	20700060

VACUU-SELECT configurations for control in fine vacuum range

VACUU-SELECT fine vacuum control packages provide an easy-to-use vacuum control option for work between 1 mbar and 10^{-3} mbar, such as freeze drying, Schlenk lines, and molecular distillation. Predefined applications can be adjusted easily, and new applications can be created in a matter of seconds, to meet your precise needs.

Each fine vacuum control package includes the VACUU-SELECT controller, fine vacuum sensor, solenoid control valve, and connections. Choose between a Pirani-Sensor or - with extended measuring range - the VACUU-VIEW extended. All valves and sensors use VACUU-BUS communication. For use with smaller fine vacuum pumps like the RE/RZ 2.5 and RE/RZ 6 rotary vane pumps, or RC 6 chemistry-HYBRID pump, choose a KF DN 16 fine vacuum control package. For larger fine vacuum pumps like the RE/RZ 9 rotary vane pump, use a KF DN 25 fine vacuum control package.




pressure sensor	control range	measuring principle	vacuum connector + Ord. No.	
VSP 3000	5 - 10^{-3} mbar	ceramic coated Pirani-Sensor	KF DN 16 / HN DN 10:	20700101
			KF DN 25 / HN DN 15:	20700111
VACUU-VIEW extended	atm. - 10^{-3} mbar	ceramic coated Pirani-Sensor + capacitance sensor	KF DN 16 / HN DN 10:	20700100
			KF DN 25 / HN DN 15:	20700110

KF = small flange HN = hose nozzle

VACUU-SELECT chemistry pumping units

These chemistry pumping units use chemical resistant solenoid valves to maintain process control. Once the first boiling point is detected by the controller, it maintains the pressure at this level. Pumping units can be configured to run one or two independent processes at the same time.

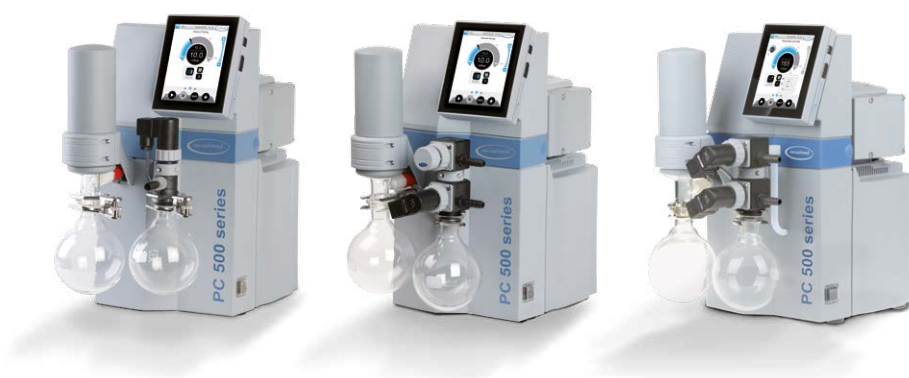

Pumping units with one electronically controlled port



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 510 select	7 mbar	2 m ³ /h	CEE: 20733150	CH: 20733151	UK: 20733152
			US: 20733153	CN: 20733156	IN: 20733157
PC 610 select	1.5 mbar	3.4 m ³ /h	CEE: 20737150	CH: 20737151	UK: 20737152
			US: 20737153	CN: 20737156	IN: 20737157

Pumping units with two vacuum ports (One electronically controlled port and one manual flow control)



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 511 select	7 mbar	2 m ³ /h	CEE: 20733250	CH: 20733251	UK: 20733252
			US: 20733253	CN: 20733256	IN: 20733257
PC 611 select	1.5 mbar	3.4 m ³ /h	CEE: 20737250	CH: 20737251	UK: 20737252
			US: 20737253	CN: 20737256	IN: 20737257

Pumping units with two electronically controlled ports



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 520 select	7 mbar	2 m ³ /h	CEE: 20733350	CH: 20733351	UK: 20733352
			US: 20733353	CN: 20733356	IN: 20733357
PC 620 select	1.5 mbar	3.4 m ³ /h	CEE: 20737350	CH: 20737351	UK: 20737352
			US: 20737353	CN: 20737356	IN: 20737357



VARIO® CONTROLLED PUMP MODELS

VARIO select chemistry diaphragm pumps

VACUU·SELECT simplifies your work and allows VARIO select chemistry diaphragm pumps to deliver unparalleled precision. The pumps run whisper quiet and respond to demand in real time by matching pumping speed to process conditions. This minimizes power consumption and maintenance. For solvent evaporation, vapor pressures are automatically detected and continuously adapted throughout the process.

The intuitive VACUU·SELECT controller includes predefined applications for all common vacuum processes. Application parameters can be adjusted quickly, and new applications can be created in seconds to suit your needs. VACUUBRAND VARIO® technology - frees you up for other tasks.



product	ultimate vacuum	max. flowrate	Ord. No.		
MZ 2C VARIO select	7 mbar	2.8 m ³ /h	CEE: 20732450	CH: 20732451	US: 20732453
MD 4C VARIO select	1.5 mbar	4.6 m ³ /h	CEE: 20736550 CN: 20736556	CH: 20736551 IN: 20736557	US: 20736553



product	ultimate vacuum	max. flowrate	Ord. No.		
ME 16C VARIO select	70 mbar	20 m ³ /h	CEE: 20741750	US: 20741753	
MD 12C VARIO select	1.5 mbar	15 m ³ /h	CEE: 20743750 CN: 20743756	UK: 20743752	US: 20743753
MV 10C VARIO select	0.6 mbar	13 m ³ /h	CEE: 20744750 CN: 20744756	UK: 20744752	US: 20744753

VARIO select chemistry pumping units

VARIO pumping units offer same benefits as VARIO diaphragm pumps. Additionally, all pumping units include an inlet catchpot (separator), an outlet catchpot, and a solvent condenser. The inlet catchpot collects particulate and fluid matter from the vacuum line. The outlet catchpot and solvent condenser combine to prevent solvent vapor emissions into the lab, allowing for nearly full recovery of solvents.



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 3001 VARIO select	2 mbar	2 m ³ /h	CEE: 20700200	CH: 20700201	UK: 20700202
			US: 20700203	CN: 20700206	IN: 20700207



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 3002 VARIO select	7 mbar	2.8 m ³ /h	CEE: 20733550	CH: 20733551	UK: 20733552
PC 3004 VARIO select	1.5 mbar	4.6 m ³ /h	US: 20733553	CN: 20733556	IN: 20733557
			CEE: 20737550	CH: 20737551	UK: 20737552
PC 3003 VARIO select	0.6 mbar	2.8 m ³ /h	US: 20737553	CN: 20737556	IN: 20737557
			CEE: 20738450	CH: 20738451	UK: 20738452
			US: 20738453	CN: 20738456	IN: 20738457



product	ultimate vacuum	max. flowrate	Ord. No.		
PC 3016 VARIO select	70 mbar	20 m ³ /h	CEE: 20741850	US: 20741853	
PC 3012 VARIO select	1.5 mbar	15 m ³ /h	CEE: 20743850	CH: 20743851	UK: 20743852
			US: 20743853	CN: 20743856	IN: 20743857
PC 3010 VARIO select	0.6 mbar	13 m ³ /h	CEE: 20744850	CH: 20744851	UK: 20744852
			US: 20744853	CN: 20744856	IN: 20744857



Email: enquiries@asynt.com **Tel:** +44 (0)1638 781 709 **Web:** www.asynt.com